

Cleaning of 25/96-well plates

1. Rinse wells with ddH₂O several times by pipetting ddH₂O into each well. Pipette up and down for each well, then shake water out by inverting the plate upside-down.
2. Repeat 3 -4 times.
3. Pipette 100 EtOH into each well and allow to sit for 5 minutes.
4. Shake out EtOH by inverting the plate upside-down and shaking.
5. Pipette 70% EtOH into each well, and allow to sit for 10 min.
6. Shake out EtOH by inverting the plate upside-down and shaking.
7. Under a sterile hood, rinse wells with ddH₂O, shake out water and allow to air dry.